IRIS Salmonella®

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(x) g of sample in 9 (x) mL of *Salmonella* Enrichment medium ②

+

Add selective IRIS *Salmonella*® supplement

Green coloration

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
Plate 10 µL on IRIS *Salmonella*® Agar

D+1

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D+1

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Reading and confirmation

Non characteristic colonies :
ABSENCE of *Salmonella*

Characteristic magenta colonies:
PRESUMPTIVE PRESENCE of
Salmonella

Latex coagulation test :
CONFIRM' *Salmonella*

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① SIRIS *Salmonella* method is also validated for the detection of *Salmonellae* in samples from 50 g to 375 g in milk powder (including infantile formula with and without probiotics) and from 50 g to 125 g for flours and croquettes in feedstuffs.

② For the samples greater than 25g, preheat the *Salmonella* Enrichment medium at 41.5° C.

Salmonella Enrichment and IRIS *Salmonella*® Agar can be stored up to 3 days at 2-8 ° C before plating or reading (except the enrichment for the sample greater than 25 g in feedstuffs).

Salmonella Enrichment can be used for the preparation of dilutions.

IRIS *Salmonella*® Agar may be used as the second medium according to NF EN ISO 6579.

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Salmonella Enrichment

BK194HA/GC – 500 g bottle/5 kg drum
BM13608 – 10 vials of 225 mL
BM13708 – 3 flexible bags 3 L
BM14408 – 2 flexible bags 5 L

Salmonella Enrichment + Tween®80

BM16308 – 3 flexible bags 3 L
BM19808 – 2 flexible bags 5 L
BM21608 – 10 vials of 225 mL

Salmonella Enrichment double-strength buffered

BM20008 – 2 flexible bags 5 L
BM20108 – 10 vials of 225 mL
BK225HA/GC – 500 g bottle/5 kg drum

IRIS *Salmonella*® supplement

BS07808 – 10 vials of 50 mL
BS07708 – 120 tablets qsp 225 mL
BS09308 – 120 tablets qsp 90 mL

IRIS *Salmonella*® Agar

BM16008 – 20 Petri dishes (Ø90mm)
BM16108 – 120 Petri dishes (Ø90mm)
BK212HA – 500 g bottle

CONFIRM' *Salmonella*

BT01108 – Latex agglutination kit
(50 tests)